

Hydraulic Arm Instruction



Attention:

Different models of stands may be received according to your purchase order, so please use the same instructions. If there is any question, please call the manufacturer or local dealer.

Step.1

Open the stand box, and check the packed contents.

The parts list

1. Clamp
2. Arm
3. Tight Screw
4. Focusing Scope Holder



Step 2. Put Clamp 1 on the side of the table and tighten the clamp.

Step 3. Insert Screw 3 inside the hole of the clamp.

Step 4 Put Arm 2 on Screw 3, and fasten the screw from the top, and cover the hole with the cap.



Step 5. The clamp has an adjustable screw on its side. Fasten the screw to tighten the arm.

Step 6. Put Focusing Scope Holder 4 in front of the Arm. Slide down the holder and fasten the two top screws, then slide up the holder and fasten the two bottom screws.



Step 7. An adjustable screw on the end of the Arm is for adjusting the tightness or looseness depending on the weight to be put on the front part of the scope holder.

Attention: When lower level position is required for the arm, press the arm down to the position, and pull up the e-arm, push down the end of the arm and keep the focusing holder at level position.



Howard Electronic Instruments, Inc. Address: 974 SE Pioneer Road. El Dorado, Kansas 67042

Tel: 1-800-394-1984 Fax: 316-321-2803 Website: www.heiscope.com Email: sales@heiscope.com